Nota:

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Highlights

- Results in 3-5 minutes
- Simple protocol

Contents of Kit

- 50 QuickTox Strips packed in a moisture-resistant canister
- DB2 Buffer
- 100 pipette tips
- 50 reaction vials
- MMBC

Items Not Provided:

- Orbital/rotary shaker
- Plastic sample cups w/lids*
- 100 μ L fixed-volume pipette* (or adjustable pipette)
- 50% laboratory grade ethanol
- 20 mesh screen
- Graduated cylinder*
- Approved coffee filter*
- Timer

* Available as accessories – see list on Page 6

Correct 20 mesh grind for corn

Measure Solvent

Catalog Number AQ 201 BGE

Part# 13298

Intended Use

This QuickTox Kit for Aflatoxin is designed to quickly extract and screen corn for the presence of total aflatoxins utilizing a solvent-based extraction. The QuickTox Kit is designed to provide a qualitative screen for aflatoxin residues at a 20 ppb cut off level in corn grain.

How the Test Works

A composite sample is first collected, ground, and extracted to solubilize any aflatoxin present. The extract is further diluted into Buffer before being run on the QuickTox test strip.

Each QuickTox Strip has an absorbent pad at each end. The protective tape with the arrow indicates which end of the strip to insert into the reaction vial. The sample extract travels up the membrane strip and is absorbed into the larger pad at the top of the strip. The portion of the strip between the protective tape and the absorbent pad at the top of the strip is used to view the reactions as described under “Interpreting the Results.”

Sample Preparation

Step 1: Determine Number and Size of Sub-samples

- Collect a composite corn sample according to your own sampling plan or USDA/GIPSA guidelines. Consult USDA/FGIS (formerly GIPSA) reference documents such as http://www.gipsa.usda.gov/fgis/handbook/BK1/BookI_2015-09-18.pdf to help design a plan that fits your needs.
- Grind samples using a mill or grinder which provides a sample such that $\geq 95\%$ passes through a 20 mesh sieve. Mix ground material thoroughly before sub-sampling.

Step 2: Extract corn sample

Important: Ensure that all samples, ethanol, buffer, and test strips are at room temperature before use

- Weigh 10-50 grams of milled sample into a disposable sample cup with lid (or other suitable vessel) and add four volumes of 50% ethanol (4 mL per gram of sample, i.e. 10 grams, add 40 mL).
- 50% Ethanol Preparation: For 100 mL, measure 50 mL 100% ethanol [reagent grade or better], pour into suitable container with cap. Add 50 mL deionized or distilled water. Cap tightly and shake to mix. Use care when uncapping. Ethanol can also be purchased at 50% concentration.
- Cap tightly and shake for 30 seconds on a mechanical shaker operated at the highest speed. Alternately, samples may be shaken vigorously by hand for 1 minute.
- Pour sample through an approved coffee filter (ACC-083) into a clean vessel. Once enough extract has passed through (no longer than 2 minutes), move filter aside to access the extract. Alternatively, allow sample to settle in extraction cup; the top (yellowish) layer containing the aflatoxin residues will be used in testing.

Step 3: Dilute corn extract with DB2 Buffer

- Using a 100 μ L fixed pipette and a new tip, transfer 100 μ L of DB2 Buffer into the reaction vial.
- Discard tip, and using a new tip, add 100 μ L of clarified extract to reaction vial.
- Mix buffer and sample extract thoroughly by pipetting up and down a few times to homogenize well.

How to Run the QuickTox Strip Test

- Allow refrigerated canisters to come to room temperature before opening. Remove the QuickTox Strips to be used. Avoid bending the strips. Reseal the canister immediately.
- Place the strip into the reaction vial containing the diluted sample extract. The arrow tape on the end of the strip should point into the reaction vial.

Shake Mechanically...

...or by Hand

Filter Extract

*Add Buffer to Vial First –
Then Add Extract; Mix Well*

Place Strip in Vial

- The sample extract will travel up the strip (flow may not be visible immediately—this is expected and normal).
- Allow the strip to develop for 5 minutes before making final assay interpretations. Negative sample results may become obvious much more quickly (2-3 minutes). Development of a Control Line within 5 minutes indicates that the strip has functioned properly. Any strip that does not develop a Control Line should be discarded. A second preparation of the extract (using a fresh 1:1 dilution) should be made and tested using another strip.
- To retain the strip, cut off and discard the bottom section of the strip covered by the arrow tape.

Interpreting the Results

Visual Interpretation

Negative Results – A sample containing aflatoxin residues of less than 10 ppb will develop 2 distinct lines in the test area. A negative test result can be interpreted as soon as a Test Line develops, generally within 2-3 minutes.

Positive Results – The QuickTox Kit for Aflatoxin Free is designed to screen for aflatoxin at levels of approximately 20 ppb or higher in corn grain. A sample containing aflatoxin residues of 20 ppb or higher will develop 1 distinct line, the Control Line. The absence of a Test Line should be interpreted as positive for aflatoxin residues. Allow the strip to develop for the full 5 minutes before concluding that the sample has tested positive for aflatoxin. Some samples containing slightly less than 20 ppb may also provide a positive result. You may wish to confirm positive results with a quantitative method to determine the precise level of contamination.

QuickScan Interpretation

Important Notes:

- The lot-specific Multi-Matrix Barcode Card (MMBC) must be scanned into the system prior to testing.
- Reading the strip on the QuickScan is required for use of the TotalHub Data Management Platform. For more information on TotalHub, contact us at <https://www.envirologix.com/contact/>.
- Sample information may be inputted before or after reading the strip.

Reading the Test:

- Place strip in the QuickScan carrier and slide carrier in. Click “Read Test” on Main Menu. Aflatoxin results are reported as “positive” or “negative”. Negative results indicate the sample contains less than 20 ppb total aflatoxins.
- Detailed instructions for use of the QuickScan System are supplied with each unit and can also be found at www.envirologix.com/support.

Kit Storage

This QuickTox Kit should be stored refrigerated. Note the shelf life on the kit box. The kit may be used in field applications; however, prolonged exposure to high temperatures may adversely affect the test results. Do not open the desiccated canister until ready to use the strips.

Cross-Reactivity

The following mycotoxins have been tested with this kit using the protocols specified herein with no false positive results.

- DON (deoxynivalenol)
- Fumonisin B₁
- Ochratoxin A
- T-2/HT-2
- Zearalenone

Available Accessories

Graduated Cylinder.. **ACC-068**
(100 mL)

MiniPet 100 µL.....**ACC-041**

5 oz. Sample Cups with
Lids..... **20-0047**
*Case of 500, for samples
up to 30g*

Coffee Filters..... **ACC-083**
(100)

Precautions & Notes

SAFETY

- Disposal of aflatoxin-contaminated materials. Follow your facility's safety procedures for disposal of samples and extracts potentially containing or known to contain aflatoxin.
- Ethanol is flammable and toxic. Avoid inhaling vapors or contact with the skin, eyes, or clothing. Wear personal protective equipment including safety glasses, nitrile gloves (not latex), a vapor mask and a lab coat when handling. Keep containers tightly closed and away from heat, sparks and open flame. Observe any applicable regulations when disposing of samples and kit reagents.

GENERAL

- This kit is designed to screen for presence or absence only, and is not designed to be quantitative.
- This product is currently not applicable for use in testing any other crops.
- As with all screening tests, it is recommended that results be confirmed by an alternate method when necessary.
- The assay has been optimized for use with the protocol provided in the kit. Deviation from this protocol may invalidate the results of the test.
- The results generated through the proper use of this diagnostic tool reflect the condition of the working sample directly tested. Extrapolation as to the condition of the originating lot, from which the working sample was derived, should be based on sound sampling procedures and statistical calculations which address random sampling effects, non-random seed lot sampling effects and assay system uncertainty. A negative result obtained when properly testing the working sample does not necessarily mean the originating lot is entirely negative for the analyte or protein in question.
- A negative result may safely be interpreted in as little as 2-3 minutes after beginning the test. It is not safe, however, to interpret positive results prior to 5 minutes.
- Protect all components from hot or cold extremes of temperature when not in use. Do not leave in direct sunlight or in vehicle.

LIMITED WARRANTY

EnviroLogix Inc. (“EnviroLogix”) warrants the products sold hereunder (“the Products”) against defects in materials and workmanship when used in accordance with the applicable instructions for a period not to extend beyond a product’s printed expiration date. If the Products do not conform to this Limited Warranty and the customer notifies EnviroLogix in writing of such defects during the warranty period, including an offer by the customer to return the Products to EnviroLogix for evaluation, EnviroLogix will repair or replace, at its option, any product or part thereof that proves defective in materials or workmanship within the warranty period.

ENVIROLOGIX MAKES NO OTHER WARRANTIES, EXPRESS OR IMPLIED, INCLUDING BUT NOT LIMITED TO ANY IMPLIED WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE. The warranty provided herein and the data, specifications and descriptions of EnviroLogix products appearing in EnviroLogix published catalogues and product literature are EnviroLogix’ sole representations concerning the Products and warranty. No other statements or representations, written or oral, by EnviroLogix’ employees, agents or representatives, except written statements signed by a duly authorized officer of EnviroLogix Inc., are authorized; they should not be relied upon by the customer and are not a part of the contract of sale or of this warranty.

EnviroLogix does not warrant against damages or defects arising in shipping or handling, or out of accident or improper or abnormal use of the Products; against defects in products or components not manufactured by EnviroLogix, or against damages resulting from such non-EnviroLogix made products or components. EnviroLogix passes on to customer the warranty it received (if any) from the maker thereof of such non-EnviroLogix made products or components. This warranty also does not apply to Products to which changes or modifications have been made or attempted by persons other than pursuant to written authorization by EnviroLogix.

THIS WARRANTY IS EXCLUSIVE. The sole and exclusive obligation of EnviroLogix shall be to repair or replace the defective Products in the manner and for the period provided above. EnviroLogix shall not have any other obligation with respect to the Products or any part thereof, whether based on contract, tort, strict liability or otherwise. Under no circumstances, whether based on this Limited Warranty or otherwise, shall EnviroLogix be liable for incidental, special, or consequential damages.

This Limited Warranty states the entire obligation of EnviroLogix with respect to the Products. If any part of this Limited Warranty is determined to be void or illegal, the remainder shall remain in full force and effect.

License

EnviroLogix has developed this kit using proprietary reagents.

EnviroLogix, the EnviroLogix logo, and QuickTox are trademarks of EnviroLogix Inc.

For Technical Support Contact Us At:

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EnviroLogix, o logotipo EnviroLogix e QuickTox são marcas comerciais da EnviroLogix Inc.